User manual
Benutzerhandbuch
Manual de usuario
Gebruikershandleiding

Microscopes

Article codes | Artikelnummern | Códigos artículos | Artikelnummers:
75895, 75925, 75938, 75939 & 75940

WWW.BMSMICROSCOPES.COM


BMS_basic-microscopes_series_GB-DE-ES-NL_manual_v1.PDF9001Basic

BMS 036 Basic
BMS 036 rechargeable
BMS 037 LED
BMS 037 LED Pro Mono
BMS 037 LED Pro Bino

GB

Thank you for choosing BMS Microscopes.

This basic compound microscope is an excellent unit for starting hobbyists. The enhanced LED illumination combines the advantages of an incident light type with a transmission type to a handy combined microscope. The transmitted and incident light are separately operable. This microscope is developed according to the latest technology on optical and mechanical details. It is recommended though that an adult is supervising at all time.

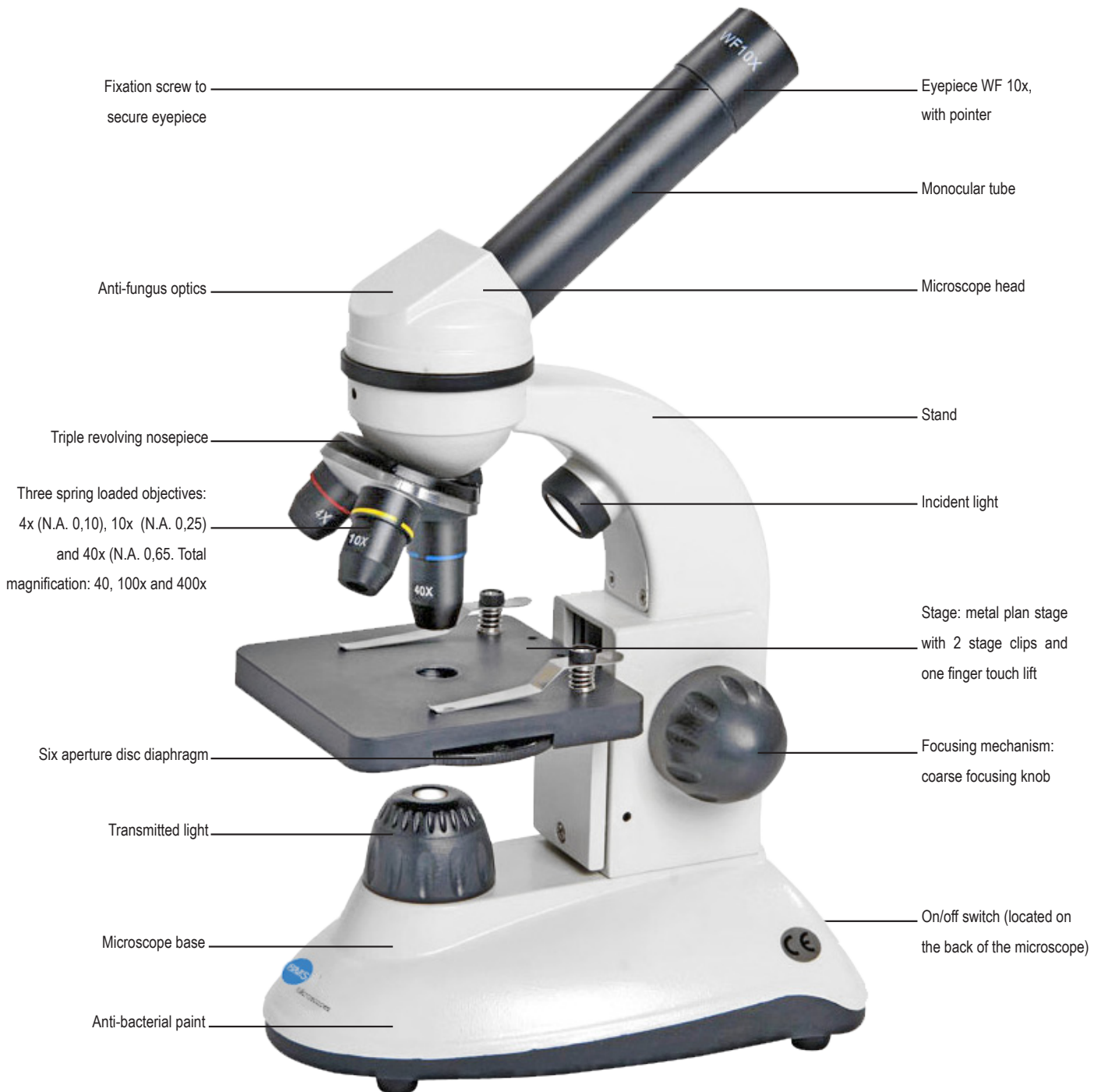Including: dust cover, manual and batteries (3x AA batteries)

Package dimensions (Width x Height x Depth): 21 x 35 x 16 cm

Gross weight: 1,1 kg

Warranty: 5 years

Article code: 75895

This basic model works with LED battery operated light source and is ideal for taking with you on the go. This very handy compound microscope is ideal for beginners! It allows you to study structures and textures of botanical and zoological nature at cell level. This microscope is rechargeable. It is developed according the latest technology on optical and mechanical details. It is recommended though that an adult is supervising at all time.

Including: dust cover, manual, 3 rechargeable AA batteries and charger

Package dimensions (Width x Height x Depth): 21 x 34 x 16 cm

Gross weigh: 1,8 kg

Warranty: 5 years

Article code: 75925

The new BMS 037 LED (Pro) is an ideal starter microscope for basic level education and hobbyists. The stability of the microscope is optimized by the robust stand and construction. Because of its cleverly placed handle in the stand, the BMS 037 LED (Pro) will move effortlessly wherever you need it to be. This microscope is developed according to the latest technology on optical and mechanical details. The enhanced LED illumination combines the advantages of an incident light type with a transmission type to a handy combined microscope. The transmitted and incident light are separately operable and are adjustable by a dimmer. This light source is characterized by low maintenance and low power consumption. This LED source has an average life span of 50,000 hours. Beside that, this LED illumination system lacks heat development and is therefore safe to use by children. It is however always recommended that an adult is supervising at all time.

Including: dust cover, manual, 3 rechargeable AA batteries and charger with charge indicator

Package dimensions (Width x Height x Depth): 28 x 38 x 20 cm

Gross weigh: 2,2 kg

Warranty: 5 years


Article code: 75938

BMS 037 Pro MONO
Article code: 75939

BMS 037 Pro BINO
Article code: 75940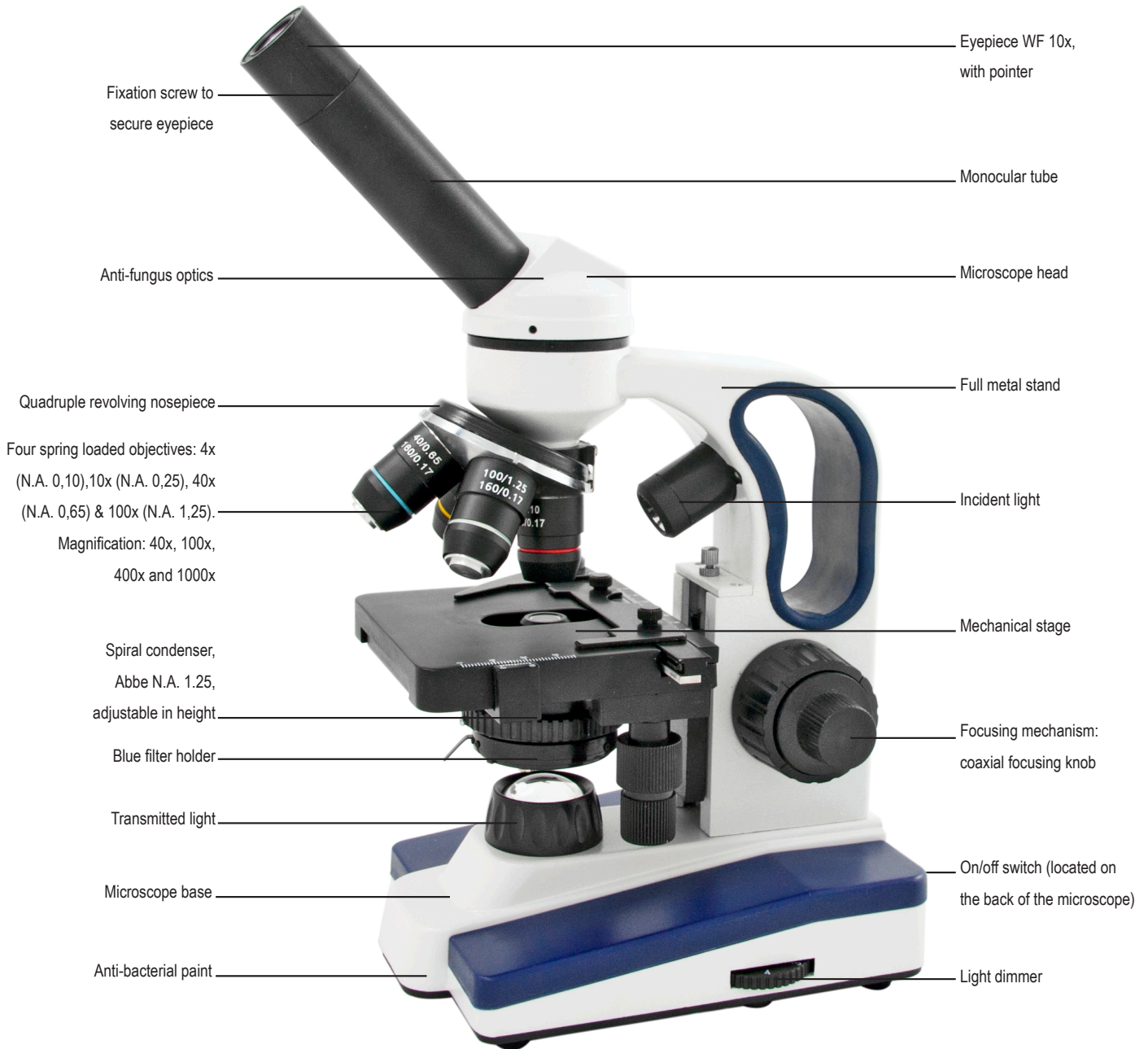

Including: dust cover, manual, 3 rechargeable AA batteries and charger with charge indicator

Package dimensions (Width x Height x Depth): 28 x 38 x 20 cm

Gross weigh: 2,6 kg (mono) and 3 kg (bino)

Warranty: 5 years

GETTING STARTED

The objectives have already been placed in the microscope. They are parfocal and parcentered. Unless necessary, do not remove the objectives. Remove the microscope carefully from the package. Open the battery holder on the bottom of the microscope. Place the batteries in the microscope. Starting with the -side and then the +side. Close the battery holder. The microscope is now ready for use.

POWER OPTIONS

1. Main power adapter.
2. Rechargeable batteries (AA, 1,5V), only applicable for rechargeable microscopes. It can be ordered as a pack of three (item no. 29554). When this battery pack is placed, it will be charged automatically when the microscope is connected to the mains and the power switch of the microscope is switched on. When the LED indicator, located on the back of the microscope, is green, the battery pack is fully charged. The microscope can now be used, disconnected from the mains, for 10 hours. When the LED indicator is red, the batteries need to be charged again. Charging time is about 3 hours.
3. Non rechargeable batteries (AA 1,5V, 3 pieces). The microscope will also work with these batteries. **WARNING:** do not connect the microscope to the main electricity when these batteries are placed. This will cause severe damage to the microscope.

SAFETY INFORMATION

- Always disconnect mains plug from power point before maintenance.
- Work in a clean and proper environment.
- Think of the safety of yourself and those working around you.
- Clean all parts coming in contact with the observed specimen to prevent infection.
- Place the microscope on a sturdy, level table or working surface.
- Always use original accessories.

CAUTION

The delicate components may get damaged if the microscope is not handled with care. Use the microscope in a clean environment. Do not use it when it is subject to direct sunlight, high temperatures, humidity or dust. Carry the microscope with care, do not hold it by the stage. Remove specimen before transport.

CARE & MAINTENANCE

- Please use our special microscope maintenance kit (item no. 76026).
- Always keep the optical parts clean. Please use the dust cover to protect it.
- Avoid contamination of the objectives.
- In case there is dirt on the lens surface, use distilled water to moist and soft cloth to clean the lens surface. Be very careful and make sure to avoid scratches on a lens surface. Cleaning can also be done by using an alcohol solution or an optical cleaning solution. Ventilate the room. Do not soak the lens in (cleaning) liquid.
- To clean the eyepiece- and objective-lens surface, move the cloth in a spiral motion from the center to the rim.
- Never make a zig-zag movement since this will only spread the dirt.
- To clean painted parts, use also a slightly moistened soft cloth.
- Never use an organic solvent. Use a mild detergent, like soft hand soap.
- Do not remove the objectives from the nosepiece. The objectives were placed in a particular sequence and are adjusted in order to have a parcentered optical system.
- Never dismantle the microscope. This can only be done by factory and/or dealer authorized staff.

SPECIFICATIONS

For more specifications, please view the product page on the website: www.bmsmicroscopes.com

It is within our company policy to develop continuously. BMS Microscopes b.v. reserves the right to change design and specifications without further notice.

BMS Microscopes b.v.

Essebaan 50

2908 LK Capelle aan den IJssel


The Netherlands

Tel.: +31(0)10 458 42 22

E-mail: sales@bmsmicroscopes.com

Environment protection first!

Your appliance contains valuable materials which can be recovered or recycled. Leave it at a local civic waste collection point.

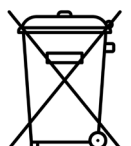General troubleshooting tables

If difficulties should be encountered in the course of operation, and no major instrument malfunctions can initially be detected, please recheck the symptoms, referring to the tables provided below, before contacting your service representative.

OPTICAL SYMPTOMS	CAUSES	COUNTERMEASURES
Darkness at the periphery, or uneven viewfield brightness	Revolving nosepiece not in click-stop position (objective not centered in optical path)	Revolve to click-stop position (swing the objective correctly into the optical path)
	Field diaphragm not centered	Centering
	Field diaphragm closed too far	Open properly
	Dirt or dust on the lens(condenser, objective, eyepiece etc)	Cleaning
Dirt or dust in the viewfield	Dirt or dust on the lens (condenser, objective, eyepiece etc)	Cleaning
	Dirt or dust on the slide	Cleaning
	Condenser position too low	Correct positioning
Poor image quality (low resolution or contrast)	No coverglass attached to the slide	Attach coverglass
	Coverglass too thick or thin	Use coverglass of specified thickness (0.17mm)
	Slide upside-down	Turn over the slide
	Immersion oil on dry system objective (especially 40X)	Cleaning
	Air bubbles in immersion oil	Remove bubbles
	Condenser aperture and field diaphragm opened too far	Close properly
	Dirt or dust on the entrance lens	Cleaning
	Condenser aperture closed too far	Open properly
	Condenser position too low	Raise to the position where the diaphragm image is in focus
Image dark on one side	Revolving nosepiece not in click-stop position	Revolve to click-stop position
	Floating specimen	Fasten securely
Image shifts during focus	Specimen rises from stage surface	Place it stable
	Revolving nosepiece not in click-stop position	Revolve to click-stop position
Image tinged yellow	Daylight filter not used	Use daylight filter
Insufficient illumination brightness	Condenser aperture too small	Readjust aperture
	Condenser position too low	Correct positioning
	Dirt or dust on the lens (condenser, objective, eyepiece etc)	Cleaning

MECHANICAL SYMPTOMS	CAUSES	COUNTERMEASURES
Specimen image unsmooth	Mechanical stage not securely fastened	Tighten all fasteners
Image not focusable with high power objectives	Slide upside-down	Turn slide over
	Coverglass too thick	Use coverglass of specified thickness (0.17mm)
Binocular image not integrated	Interpupillary distance not correctly adjusted	Adjustment
Excessive eye fatigue	Incorrect diopter adjustment	Correct adjustment
	Inadequate brightness or illumination	Adjust brightness with control dial

ELECTRICAL SYMPTOMS	CAUSES	COUNTERMEASURES
Lamp does not light when switched ON	No electrical power	Check power cord connection
	LED Lamp not inserted	Insert correctly
	LED failure	Replacement: if the same symptom occurs after replacing the lamp, contact your dealer
Insufficient illumination brightness	Unspecified lamp bulb used	Replace with specified lamp bulb
	Voltage too low	Increase brightness with control dial
Unstable lamp brightness	LED Lamp about to fail	Replacement
	LED Lamp not correctly inserted into socket	Check for positive connection

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